

WESTSAVE Up™ (Western Blotting Substrate)

For research use only


Description

AbFrontier WESTSAVE Up™ (Western Blotting Substrate) is enhanced luminol-based chemiluminescent substrate for the non-radioactive detection of Horseradish Peroxidase(HRP) labeled antibodies.

Contents

LF-QC0101	WESTSAVE Up™ (Western Blotting Substrate)	
	Detection reagent A (Luminol Enhancer solution)	200 ml
	Detection reagent B (Peroxide solution)	1 ml
LF-QC0102	WESTSAVE Up™ (Western Blotting Substrate, double pack)	
	Detection reagent A (Luminol Enhancer solution) X 2 ea	400 ml
	Detection reagent B (Peroxide solution) X 2 ea	2 ml

Basic Principle

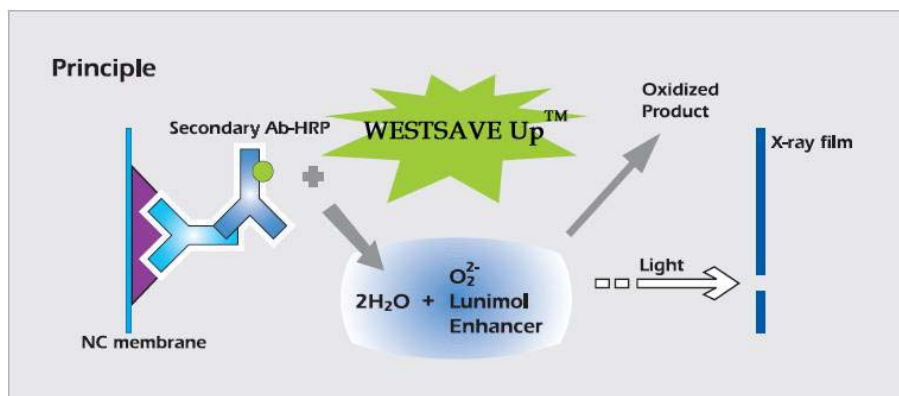

Short protocol

- Mix 5 ml of **Detection Reagent A** and 10 μl of **Detection Reagent B (500:1)**.
 - If you want to see the strong signal, you can add more 1.5~2 folds Detection Reagent B.
 - Make this mixture right before you intend to use the solution.
- Incubate blot with mixed solution for 10 seconds by placing in a tray and drain excess reagent.
- Wrap in clear plastic sheet and then expose blot to X-ray film.

Storage: Stable for 6 months at 4°C from date of shipment.

Caution: Do not use any solution with sodium azide. It will destroy the HRP activity.
Wash buffer(pH8.0) is recommended.

Abfrontier. Co. Ltd.

Tel : +82-2-6933-6789

E-mail: orders@abfrontier.com

Address: 1 Fl., Science Building C, Ewha University Daehyun-dong, Seodaemoon-gu, Seoul, 120-125, Korea

Fax: +82-2-6933-6780

Website: <http://www.abfrontier.com>